

Mut Express MultiS Fast Mutagenesis Kit V2

C215



Nanjing Vazyme Biotech Co.,Ltd.

Tel: +86 25-83772625

Email: info.biotech@vazyme.com

Web: www.vazyme.com

Loc: Red Maple Hi-tech Industry Park, Nanjing, PRC

Follow Us



Instruction for Use

Version 22.1

Contents

01/Product Description	02
02/Components	02
03/Storage	02
04/Applications	02
05/Self-prepared Materials	02
06/Notes	03
07/Mechanism & Workflow	03
08/Experiment Process	04
08-1/Primer Design for Multi-base Site-directed Mutagenesis	04
08-2/Amplification of Target Plasmid	04
08-3/ <i>Dpn</i> I Digestion of Amplification Products	05
08-4/ <i>Dpn</i> I-digested product Concentration Determination	06
08-5/ <i>Dpn</i> I-digested product Input	06
08-6/Recombination	07
08-7/Transformation	07
08-8/Recombinant Product Identification	07
09/FAQ & Troubleshooting	08

01/Product Description

Mut Express MultiS Fast Mutagenesis Kit V2 is a kit based on ClonExpress rapid cloning technology, which can simultaneously introduce site-directed mutagenesis to 3 - 5 discontinuous sites on the target plasmid. The kit integrates the Phanta Max Super-Fidelity DNA Polymerase amplification system and the ClonExpress rapid cloning system. With superior fidelity, Phanta Max significantly reduces the possibility of introducing novel mutations during amplification. Its excellent long fragment amplification ability is widely applicable to the amplification of any plasmid less than 20 kb. The ClonExpress rapid cloning system replaces the conventional annealing-cyclization reactions with the more efficient homologous recombination reaction. Site-directed mutagenesis using the Mut Express MultiS Fast Mutagenesis Kit V2 allows for more flexible primer design and extremely low template input, thereby facilitating complete degradation of the initially methylated template. The kit is equipped with the highly optimized recombination reaction buffer and the recombinase Exnase MultiS, which specially optimized for multi-base site-directed mutagenesis. If the amplification product is highly specific, its *Dpn* I-digested product can be used directly in the recombination reaction without DNA purification. This product is the preferred kit for DNA multiple point mutation.

02/Components

Components	C215-01 (10 rxns)	C215-02 (25 rxns)
2 × Max Buffer	1.25 ml	3 × 1.25 ml
dNTP Mix (10 mM each)	50 µl	125 µl
Phanta Max Super-Fidelity DNA Polymerase	50 µl	125 µl
<i>Dpn</i> I (10 U/µl)	50 µl	125 µl
5 × CE MultiS Buffer	40 µl	100 µl
Exnase MultiS	20 µl	50 µl

03/Storage

Store at -30 ~ -15°C and transport at ≤0°C.

▲ Please avoid repeated freezing and thawing.

04/Applications

- ◇ Mutation: 3 - 5 discontinuous mutation sites (multiple bases within 50 bp apart can be counted as 1 mutation site)
- ◇ Deletion: delete a specific sequence on the vector
 - ▲ When using this kit for site-directed mutagenesis of plasmids, use host strains without defects in methylase to amplify the original plasmid!

05/Self-prepared Materials

Original plasmids and mutant primers.

Competent cells: chemically competent cells prepared from clonal strains.

Other materials: ddH₂O, PCR tubes, PCR instrument, etc.

06/Notes

For research use only. Not for use in diagnostic procedure.

07/Mechanism & Workflow

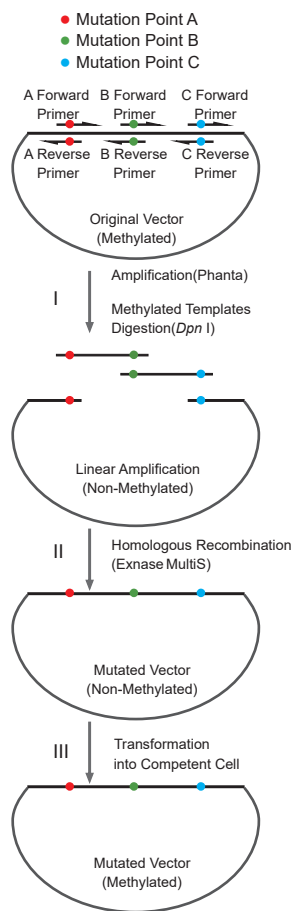


Fig 1. Flow Diagram of Discontinuous Three-base Site-directed Mutagenesis

- I. Amplification to introduce mutations: design reverse complement primers at Mutation Sites A, B and C. Use three pairs of primers for cross amplification of AF + BR to obtain fragment AB, BF + CR to obtain fragment BC and CF + AR to obtain fragment CA. Digest the amplification product using *Dpn* I.
- II. Recombination reaction: mix fragments AB, BC and CA proportionally and react at 37°C for 30 min under the catalysis of Exnase MultiS to complete the recombination reaction, achieving in vitro cyclization of linearized DNA.
- III. Transformation to competent cells: transform the recombinant product directly, and hundreds of clones will grow on the plate for positive screening later.

08/Experiment Process

08-1/Primer Design for Multi-base Site-directed Mutagenesis (Using Discontinuous Three-base Site-directed Mutagenesis as an Example)

Introduce site-directed mutagenesis into three discontinuous sites of the plasmid. It is only need to design three pairs of primers to amplify the plasmids respectively.

The primers are designed as the following: **5' - 15 - 21 bp reverse complementary region + at least 15 bp non-complementary region - 3'**

- ▲ The optimal GC content in the reverse complementary region ranges between 40% - 60%. Regions with repetitive sequences should not be selected. The optimal Tm value of the region between the target mutation site and the 3' end of the primer is >60°C. The region between the target mutation site and the 5' end of the primer should not be included in the calculation.
- ▲ The target mutation sites can be in the complementary region (point mutation should be introduced on both primers) or in the non-complementary region of either primer (point mutation is only required to be introduced on one primer). The mutation site should not be at the end of the primer.

It is recommended to use the primer design software CE Design for automatic primer generation, which can be downloaded from Vazyme's official website (<http://www.vazyme.com>). For manually design, please refer to the principle below:

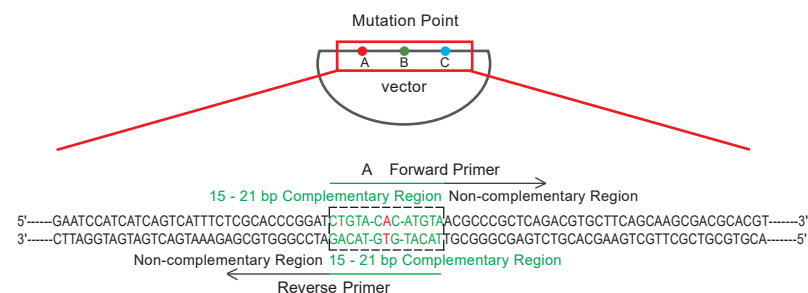


Fig 2. Primer Design for Discontinuous Multi-base Site-directed Mutagenesis

08-2/Amplification of Target Plasmid

Amplify the target plasmid using the Phanta Max Super-Fidelity DNA Polymerase. Thoroughly mix the thawed components of the reaction system by shaking. Reagents should be returned to -20°C storage immediately after use. Recommended reaction system:

Components	Volume
Template DNA ^a	Optional
2 × Max Buffer	25 μl
dNTP Mix (10 mM each) ^b	1 μl
Primer 1 (10 μM)	2 μl
Primer 2 (10 μM)	2 μl
Phanta Max Super-Fidelity DNA Polymerase ^c	1 μl
ddH ₂ O	Up to 50 μl

- Please ensure that the primers and templates do not contain uracil. And do not use dUTP.
- Use the minimum possible amount of input template if it does not affect the amplification result. It is recommended to use ≤1 ng of freshly extracted plasmids as the template.
- The recommended final concentration of the enzyme is 1 U/50 μl reaction. The optimal concentration of Phanta Max Super-Fidelity DNA Polymerase ranges between 0.5 - 2 U/50 μl, but should not exceed 2 U/50 μl, especially if the size of products >5 kb.

The recommended PCR program is as follows:

Steps	Temperature	Time	Cycles
Initial Denaturation	95°C	30 sec	1
Denaturation	95°C	15 sec	} 30 ^c
Annealing ^a	60 ~ 72°C	15 sec	
Extension ^b	72°C	30 - 60 sec/kb	
Final Extension	72°C	5 min	1

- Phanta Max Super-Fidelity DNA Polymerase promotes the efficient annealing of templates and primers. In general, the annealing temperature is set at the T_m of primers. If required, temperature gradients can be established to determine the optimal temperature for primer-template binding. A prolonged annealing time may cause diffusion of the amplification product on the gel.
- Increasing extension time can improve amplification yield.
- To prevent the introduction of non-target mutations during amplification, it is strongly recommended that the number of amplification cycles should be ≤35. If the amplification efficiency is superior, it is recommended that the number of amplification cycles should be ≤30.

After the reaction, take a small amount of amplification products for agarose gel electrophoresis detection. If the target plasmid is correctly amplified, proceed to the next step.

08-3/*Dpn* I Digestion of Amplification Products

Since the amplification product from 08-2/Amplification of Target Plasmid contains the original template plasmid, to prevent it from forming a false-positive transformant after transformation, the amplification product needs to be digested by *Dpn* I to remove the methylated template plasmid before recombination reaction. Recommended reaction system:

Components	Volume
<i>Dpn</i> I	1 μl
Amplification Product	40 - 50 μl

Please pipette up and down to mix thoroughly, then centrifuge it briefly to the bottom of the tube. Incubate at 37°C for 1 - 2 h.

- ▲ If the amplification product from step 08-2 is specific, the *Dpn* I-digested product can be used directly in the recombination reaction without purification, but the total input volume should not exceed 1/5 of the total reaction volume. If the amplified product is not specific, purify the target amplification product by gel extraction after *Dpn* I digestion.

08-4/*Dpn* I-digested product Concentration Determination

Concentration determination:

If the *Dpn* I-digested product has been purified using a high-quality gel extraction kit and no obvious non-specific bands or smears are observed in electrophoresis, its concentration can be determined using spectrophotometry-based instruments such as Onedrop. However, the concentration measured is reliable only when the A260/A280 is within the range of 1.8 - 2.0. Recommended instruments for concentration determination: Nanodrop, Onedrop, Qubit, PicoGreen.

08-5/*Dpn* I-digested product Input

The optimal amount of DNA for the recombination reaction of Exnase MultiS discontinuous multi-base site-directed mutagenesis is 0.03 pmol. The corresponding molar mass of DNA can be roughly calculated according to the following formula:

The optimal mass of *Dpn* I-digested products = [0.02 × number of fragment base pairs] ng (0.03 pmol)

For example, if fragment AB is 1 kb, fragment BC is 2 kb, and fragment CA is 5 kb, the optimal *Dpn* I-digested product input should be: 0.02 × 1,000 = 20 ng for fragment AB; 0.02 × 2,000 = 40 ng for fragment BC; 0.02 × 5,000 = 100 ng for fragment CA.

- Both excessive and insufficient DNA input amounts will reduce the cyclization efficiency. Please prepare the reaction system as strictly as possible according to the recommended amount.
- The input amount of *Dpn* I-digested products should range between 20 - 200 ng. When the optimal DNA input amount calculated using the above formula falls outside this range, select the minimum or maximum input amount within the range.
- If the *Dpn* I-digested products are used directly in the recombination reaction without purification, the total input amount should not exceed 1/5 of the reaction volume, that is 4 μl.

08-6/Recombination

The unique primer design of Mut Express (refer to 08-1/Primer Design Guidance) allows the amplified and digested products to be efficiently recombined at the target mutation region under the catalysis of Exnase MultiS to achieve in vitro cyclization of linearized DNA.

- For discontinuous multi-base site-directed mutagenesis, prepare the following reaction system on ice:

Components	Recombination Reaction	Negative Control ^b
n <i>Dpn</i> I-digested products ($n \leq 5$) ^a	$X_1 - X_n$ μ l	$X_1 - X_n$ μ l
5 \times CE MultiS Buffer	4 μ l	0 μ l
Exnase MultiS	2 μ l	0 μ l
ddH ₂ O	Up to 20 μ l	Up to 20 μ l

- X is the amount of each digestion product calculated according to the formula.
- Excessive amounts of amplification templates or incomplete *Dpn* I digestion is likely to result in a high false positive rate, so a negative control is recommended.

- Mix the reaction system thoroughly by gently pipetting (DO NOT VORTEX!), and centrifuge it briefly to collect the reaction solution to the bottom of the tube.
- Incubate at 37°C for 30 min and immediately chill the tube at 4°C or on ice.
 - ▲ It is recommended to perform the reaction on an instrument with precise temperature control such as the PCR machine. The recombination efficiency reached the highest at about 30 min of reaction. Insufficient or too long incubation time will reduce cloning efficiency.
 - ▲ The recombinant product can be stored at -20°C for one week. Thaw and transform when needed.

08-7/Transformation

- Thaw the competent cells on ice.
- Pipette 10 μ l of the recombination products to 100 μ l of competent cells, flick the tube wall to mix thoroughly (DO NOT VORTEX!), and then place the tube on ice for 30 min.
 - ▲ The volume of recombination products should be $\leq 1/10$ of the volume of competent cells.
- Heat shock at 42°C water bath for 45 sec and then immediately place on ice for 2 - 3 min.
- Add 900 μ l of SOC or LB liquid medium (without antibiotics). Then, shake at 37°C for 1 h at 200 - 250 rpm.
- Preheat the corresponding resistant LB solid medium plates in a 37°C incubator.
- Centrifuge the culture at 5,000 rpm (2,400 \times g) for 5 min, discard 900 μ l of supernatant. Then, use the remaining medium to suspend the bacteria and use a sterile bent glass rod to gently spread on the plate which contains the appropriate selection antibiotic.
- Incubate at 37°C for 12 - 16 h.

08-8/Recombinant Product Identification

After overnight incubation, if the number of clones on the plate is significantly higher than that of the negative control, select and incubate several single clones in the LB liquid medium with appropriate antibiotics overnight, and then extract the plasmids for sequencing.

09/FAQ & Troubleshooting

◇ The plasmid template cannot be amplified normally

- It is recommended to use the software CE Design for primer design. Select the appropriate module for primer design.
- Incorrect primer design: Check the primer design guidance.
- Incorrectly prepared amplification system: Repeat the experiment.
- Inappropriate amplification reaction conditions: Adjust the Mg²⁺ concentration, the amount of enzyme, and the amplification program.
- Low quality template plasmid: Long-term storage and repeated freeze-thaw cycles may cause breakage, nicking, or degradation of template plasmids. Therefore, use freshly prepared plasmids as the template.

◇ Few or no clones are formed on the plate

- Add insufficient or excessive amounts of DNA or not an appropriate ratio: Please use the amount and ratio according to specification recommended.
- Impurity of DNA inhibits the recombination reaction: The total volume of unpurified DNA should be ≤ 4 μ l (1/5 of the total volume of reaction system). It is recommended that the linearized vector and PCR product are purified by gel extraction. Then, dissolve the purified product in ddH₂O (pH 8.0).
- The low efficiency of competent cells: Make sure the transformation efficiency of competent cells is $> 10^8$ cfu/ μ g. The simple test can be performed. Transform the 0.1 ng of plasmids and take the 1/10 for spreading plates. If 1,000 clones are grown, the estimated transformation efficiency is 10^8 cfu/ μ g. The transformation volume of recombinant products should be $\leq 1/10$ of the volume of competent cells, otherwise the transformation efficiency will be reduced. Select competent cells used for cloning. Do not select competent cells used for expression.

◇ Site-directed mutagenesis is not done correctly

- Incorrect primer design: Check the primer design guidance.
- The templates used for amplification reaction are not methylated: *Dpn* I can only recognize methylated DNA. Please use the plasmids amplified from the host strain without defects in methylase as the PCR template.
- Excessive template plasmids used in the amplification reaction: For most plasmids, 1 ng of input template is sufficient for the amplification. Excessive amounts of template will lead to incomplete *Dpn* I digestion and reduce the success rate of mutagenesis.

◇ Mutations at non-target sites

- ① Template plasmids carrying unknown mutations: Check if the template plasmid sequence is correct by sequencing.
- ② Excessive amplification cycles: To prevent non-target mutations during amplification, the number of amplification cycles should be ≤ 35 . If the amplification efficiency is good, the recommended number of amplification cycles should be ≤ 30 .

◇ Special reminder!

When selecting the reverse complementary region of the primer, the region with repeated sequences should be avoided. When the GC content of this region is in the range of 40% - 60%, the efficiency of recombinant reaction will reach the maximum. If the GC content of this region is higher than 70% or lower than 30%, the efficiency of recombinant reaction will be greatly affected.